



Instrument Engineering Change Notice

NO. TEC6-2104

Sakura Finetek Japan Co. Ltd.

PRODUCT NAME : Tissue-Tek TEC6 EM DATE : December 11, 2021

PRODUCT CODES : 5104, 5108, 5112, 5116, 5128 ORIGINATOR : Yoshizawa

IECN TYPE : Part Design Change REFERENCE NUMBER : 17137-1, 17138-1

SERIAL NUMBERS AFFECTED : See Remarks ATTACHMENTS : N/A

FIELD RETROFIT REQUIRED : NO

Title: Change of Heat-shrinkable Tube in Heater Units

[REASON AND DESCRIPTION FOR CHANGE]

The heat-shrinkable tube used in heater units was changed to improve work efficiency. However, new heater units need a new fuse retainer. This makes the old and new heater units not interchangeable with each other.

Therefore, the heater unit replacement kits are available for replacement with old heater units. Attach the new fuse retainer to the new heater unit at replacement.

[PART CHANGE] (IVDP=Inventory Disposal, ICHB=Inter Changeable)

Description	Old	QTY	IVDP	New	QTY	ICHB
Fuse Retainer	A-MAL2-0202-01-A	1	NO	---	---	NO
Fuse Retainer 2	---	---		A-MAL2-0221-00-A	1	NO
Dispensing Block Heater Unit 100/115V / E:TEC6-036	A-EAL2-U036-00-A	1	NO	---	---	NO
Dispensing Block Heater Unit 100/115V / E:TEC6-056	---	---		A-EAL2-U056-00-A	1	NO
Dispensing Block Heater Unit 230V / E:TEC6-046	A-EAL2-U046-00-A	1	NO	---	---	NO
Dispensing Block Heater Unit 230V / E:TEC6-066	---	---		A-EAL2-U066-00-A	1	NO
Fuse Retainer	A-MAL2-0170-00-A	1	NO	---	---	NO
Fuse Retainer 1	---	---		A-MAL2-0171-00-A	1	NO
Forceps Holder Heater Unit 100/115V / E:TEC6-038	A-EAL2-U038-00-A	1	NO	---	---	NO
Forceps Holder Heater Unit 100/115V / E:TEC6-058	---	---		A-EAL2-U058-00-A	1	NO
Forceps Holder Heater Unit 230V / E:TEC6-048	A-EAL2-U048-00-A	1	NO	---	---	NO
Forceps Holder Heater Unit 230V / E:TEC6-068	---	---		A-EAL2-U068-00-A	1	NO

Heater Unit Replacement Kit:

Dispensing Block Heater Unit Replacement Kit (100/115V) - 1 x A-MAL2-0221-00-A fuse retainer 2 - 1 x A-EAL2-U056-00-A dispensing block heater unit 100/115V, E:TEC6-056	A-MAL2-U089-00-A
Dispensing Block Heater Unit Replacement Kit (230V) - 1 x A-MAL2-0221-00-A fuse retainer 2 - 1 x A-EAL2-U066-00-A dispensing block heater unit 230V, E:TEC6-066	A-MAL2-U090-00-A
Forceps Holder Heater Unit Replacement Kit (100/115V) - 1 x A-MAL2-0171-00-A fuse retainer 1 - 1 x A-EAL2-U058-00-A forceps holder heater unit 100/115V, E:TEC6-058	A-MAL2-U091-00-A
Forceps Holder Heater Unit Replacement Kit (230V) - 1 x A-MAL2-0171-00-A fuse retainer 1 - 1 x A-EAL2-U068-00-A forceps holder heater unit 230V, E:TEC6-068	A-MAL2-U092-00-A

[REMARKS]

Implementation:

J0: 5104 0300 and after

A1: 5108 0543 and after

E2: 5112 0374 and after

JC2: 5116 0141 and after

J2: 5128 0123 and after

[DISTRIBUTION]SFA SFJ SFE SCN **[APPROVAL]**

Name : T. Sugawara

Signature :

