



# Instrument Engineering Change Notice

NO. Prisma Plus-2001

Sakura Finetek Japan Co. Ltd.

PRODUCT NAME	: Tissue-Tek Prisma Plus	DATE	: March 18, 2020
PRODUCT CODES	: 6170, 6171, 6172, 6173, 6174, 6175, 6177	ORIGINATOR	: A. Ito, T. Iijima
IECN TYPE	: SOFTWARE CHANGE AND PART DESIGN CHANGE	REFERENCE NUMBER	: 16828-1, 16827-1, 16803, 16720
SERIAL NUMBERS AFFECTED	: See Remarks	ATTACHMENTS	: A-SK25-0900-05-A.zip
FIELD RETROFIT REQUIRED	: NO		

## Title: Software and Hardware Changes for New CI

### [ REASON AND DESCRIPTION FOR CHANGE ]


- (1) The logo displayed at the top left corner of each operation screen was deleted due to the change to the new CI logo.

The loading screen was also changed with the change of the initial program loader, as follows:

\* The initial program loader is not supplied for field service.



New loading screen (white background)

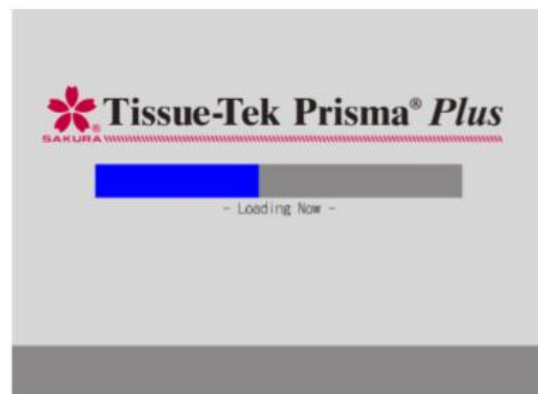

Old loading screen



New operation screen



Old operation screen

- (2) The logo on the front monitor cover was changed to the new CI logo. The cover with the old logo is supplied as a service part.  
 In addition, the monitor unit (old logo) is newly available as a service part, consisting of all the components including the control panel, CPU board, display monitor, monitor arm and monitor cover with the old logo.

**[ PART CHANGE ] ( IVDP=Inventory Disposal, ICHB=Inter Changeable )**

Description	Old	QTY	IVDP	New	QTY	ICHB
Main Control Program	A-SK25-0900-04-A	1		A-SK25-0900-05-A	1	YES
Monitor Cover F (CI)	---	---		A-MK25-0504-00-A	1	NO
Monitor Cover F	A-MK25-0500-01-A	1	NO	---	---	NO
Monitor Unit	---	---	NO	A-MK25-U012-00-A	1	NO

(Note) If the new main control program A-SK25-0900-05-A is installed to the system with an older version of the program, the loading screen is NOT updated although the operation screen is updated. This does not affect functions and performance of the system.

**[ REMARKS ]**

AD 6170 /AS 6171: K.D. (from SFA PO#202054)  
 ED: 6172 0171 and after  
 ES: 6173 0226 and after  
 JD: 6174 0250 and after  
 JS: 6175 0067 and after  
 JCS: 6177 0144 and after

**[ DISTRIBUTION ]**

SFA   
 SFJ   
 SFE

**[ APPROVAL ]**

Name : T. Sugawara

Signature :

