



Instrument Engineering Change Notice

NO. VIP6 AI-2004

Sakura Finetek Japan Co. Ltd.

PRODUCT NAME : Tissue-Tek VIP6 AI DATE : October 21, 2020
PRODUCT CODES : 6040, 6041, 6042, 6043 ORIGINATOR : Sakurai, Yokomizu
IECN TYPE : NEW PART REFERENCE NUMBER : 16984, 17130
SERIAL NUMBERS AFFECTED : See Remarks ATTACHMENTS : N/A
FIELD RETROFIT REQUIRED : NO

Title: Application of Silicon Caulking for the LCD Module

[REASON AND DESCRIPTION FOR CHANGE]

If solvents are attached to the touch panel screen, it may run down its surface down to the FPC cable terminal which may eventually cause instrument failure. To prevent the problem, silicon caulking is applied to the FPC terminal for protection from the July/August 2020 production lot.

For replacement of non-caulked LCD modules, the caulked LCD module unit is available as a maintenance part. Note that the display surface of the new caulked LCD module unit is already covered with the touch panel protection sheet (O70-626-00).

[PART CHANGE] (IVDP=Inventory Disposal, ICHB=Inter Changeable)

Description	Old	QTY	IVDP	New	QTY	ICHB
LCD Module Unit, caulked	---	---		A-MO79-U052-00-A	1	YES

[REMARKS]

Implementation:
A1: 6040 0771 and after
J0: 6041 0504 and after
E2: 6042 0773 and after
J2: 6043 0232 and after

[DISTRIBUTION]

SFA
SFJ
SFE
SCN

[APPROVAL]

Name : T. Sugawara

Signature :